

MP250M

The motor power hydraulic press

Easiest handling for frequent use

The MP250M motor power press is made to handle a large number of press actions.

You control the load by positioning a red mark inside the 63 mm pressure gauge. Although it is possible to set 2 tons as minimum load, we recommend to use die sets not smaller than 8 mm because the 63 mm pressure gauge scale is 1 ton. So the load is reproducible as long as the mark is at the same position, but it is not exact enough for very small dies. The MP250M is able to hold the load for several hours.

We integrated two additional safety features compared to the manual versions: a safety switch to prevent press actions with opened cover and an illuminated emergency stop button.

The handling is very easy: pushing the green button starts the action and it stops at the marked position. The manual release valve allows a smooth pressure decrease to avoid cracks in your pellets.



Model	MP250M
Dimensions (H min/max x W x D)	497/609 x 360 x 400 mm
Width between columns	150 mm
Min. / Max. inner height for die set	73 mm / 185 mm
Display / resolution	Analogue, 0.5 to.
Top Spindle diameter	45 mm
Max. way of spindle movement	112 mm
Cylinder diameter / max. lifting	105 mm / 22 mm
Net. weight	68 kg
Mains connection	230 V / 50 Hz / 250 W
Order Number	54MP250M



JASCO France
17 Rue des Cerisiers
91090 Lisses - France

Tél. 01 64 97 09 60 Fax: 01 85 09 90 76
email: jascofrance@jascofrance.fr